NOTICE OF APPROVAL

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Issued by statutory authority of the Minister of Industry
for:

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Chromatographe en phase gazeuse

APPLICANT

Norwest Soil Research Ltd.
7217 Roper Road
Edmonton, Alberta
T6B 3J4

REQUÉRANT

MANUFACTURER

Agilent Technologies, Inc.
2850 Centerville Road
Wilmington, Delaware
19808-1610

FABRICANT

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RATING/ CLASSEMENT

Agilent 6890A (G1540A) Gas
Chromatograph/
Chromatographe en phase gazeuse
Agilent 6890A (G1540A)

20°C to/à 27 °C

Hewlett Packard 5890-1 Series II Gas
Chromatograph
Hewlett Packard 5890-1 Série II
Chromatographe en phase gazeuse

20°C to/à 27 °C

Hewlett Packard 5890A-1 Gas
Chromatograph/Chromatographe en
phase gazeuse

20°C to/à 27 °C

NOTE: This approval applies only to meters, the design, composition, construction and performance of which are, in every material respect, identical to that described in the material submitted, and that are typified by samples submitted by the applicant for evaluation for approval in accordance with sections 13 and 14 of the Electricity and Gas Inspection Regulations. The following is a summary of the principal features only.

SUMMARY DESCRIPTION:

The Agilent 6890A (G1540A) Series or the Hewlett Packard 5890-1 Series II Gas Chromatograph analyzes the composition of natural gas. The gas chromatograph separates the following components of natural gas:

S	oxygen
S	nitrogen
S	methane
S	carbon dioxide
S	ethane
S	propane
S	iso-butane
S	n-butane
S	iso-pentane
S	n-pentane
S	hexanes
S	heptanes
S	octanes
S	nonanes
S	decanes

The individual component concentrations are used to calculate the relative density and heating value of the sample. The detection of components is performed by a TCD (thermal conductivity detector) and an FID (flame ionization detector) which provide an output signal proportional to the concentration.

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Natural gas is injected into the sampling valve and is carried by the carrier gas, helium through four columns, inside the oven. The different components move through the columns at different rates, thereby allowing identification based on the retention time of each component.

The operator interface consists of a display, status board, and keyboard on the face of the gas chromatograph as well as a user interface program (Chemstation software) running on a personal computer. The Chemstation software analyses the data and displays the chromatogram. The Chemstation software version number can be viewed by selecting the "Help" menu and then "About" in the Chemstation software program.

A Norwest Soil in house software application entitled NorLims (Norwest laboratory information management system) is used to determine heating value and relative density, using GPA 2172 (96) and GPA 2145(00) at 15°C and 101.325 kPa. The NorLims program version can be viewed under the "Help", and then the "About" menus.

The Agilent 6890A or the Hewlett Packard 5890-1 Series II gas chromatograph's firmware version can be viewed by powering down the gas chromatograph and then powering up again. The firmware version will appear on the screen during boot up.

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A Hewlett Packard model 5890A-1 analyser can be used to improve the accuracy of the gas analysis and hence the relative density and heating value calculations. Because the 6890 A or the 5890-1 Series II uses helium as a carrier gas it can not be used to detect helium so detecting helium (and hydrogen) with the 5890A-1 and then feeding that result into the NorLims software along with the output from the model 6890A (or 5890-1 Series II) enhances the analysis by making the heating value and the relative density figures more reliable. The model 5890A-1 is used only to measure these two gases and only when they are anticipated components in the gas sample under analysis. The model 5890A-1 uses only a Thermal Conductivity Detector (TCD).

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SPECIFICATIONS

Gas Chromatograph Agilent Series 6890A

Detector Type:	TCD and FID
Column 1:	6' x 1/8" - 35% DC200/500 on Chromosorb P/AW - 80/100 Back Flush Packed
Column 2:	6' x 1/8" - Haysep Q - 80/100 Packed
Column 3:	10' x 1/8" - Mole Sieve Packed 13x60/80

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MARKING REQUIREMENTS

Marking shall be in accordance with Section 3.5 and 20-3.1 of LMB-EG-08.

The model number of the HP 5890 Series II chromatograph is now 5890-1 Series II to identify it as part of the measurement system covered by this approval.

The model number of the HP 5890 chromatograph is now 5890A-1 to identify it as part of the measurement system covered by this approval.

REVISIONS**AG-0489 Rev. 1**

Revision 1 approved the Hewlett Packard model 5890A-1 chromatograph, calibrated for helium and hydrogen only, to enhance compositional resolution of the model 6890 A chromatograph.

AG-0489 Rev. 2

Revision 2 approves the Hewlett Packard model 5890-1 Series II Gas Chromatograph as it is used in the Norwest Soils Research Ltd system as an alternative to the Agilent 6890A for the primary chromatograph.

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AG-0489 Rev. 3

Revision 3 approves the NorLims analytical software application. The model number of the 5890 gas chromatograph has also been changed to 5890A-1.

EVALUATED BY**AG-0489 and Revision 3**

Judy Farwick
Complex Approvals Examiner
Tel: (613) 946-8185
Fax: (613) 952-1754

AG-0489 Revision 1 and 2

Graham Collins
Approval Examiner
Tel (613) 941-0605
Fax (613) 952-1754

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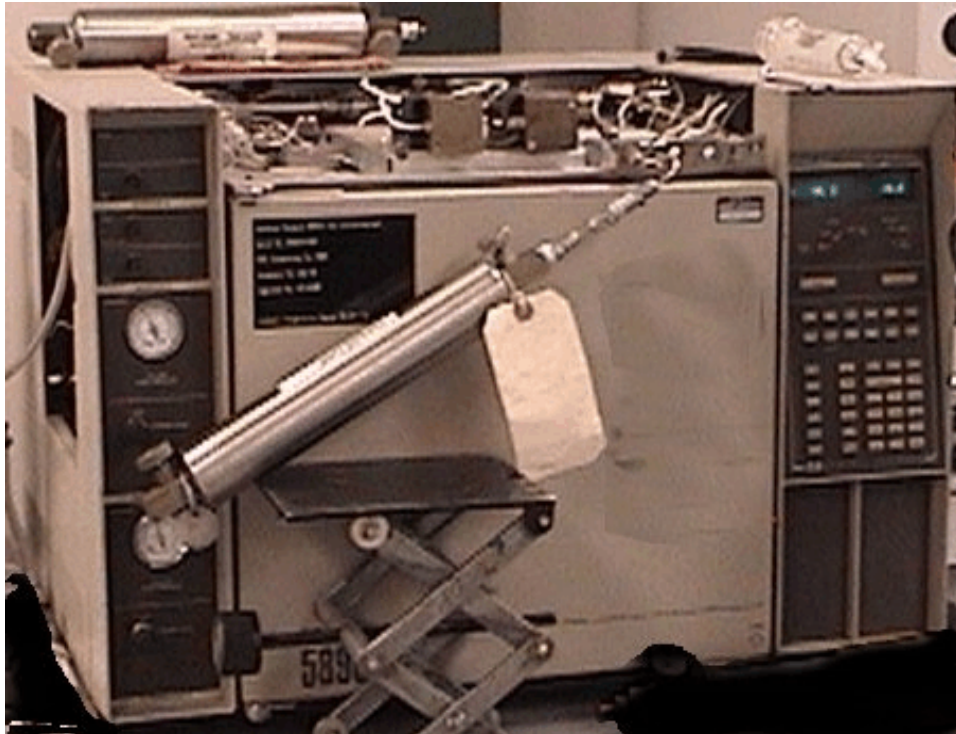
APPROVAL:

The design, composition, construction and performance of the meter type(s) identified herein have been evaluated in accordance with regulations and specifications established under the Electricity and Gas Inspection Act. Approval is hereby granted accordingly pursuant to subsection 9(4) of the Act.

The sealing, marking, installation, use and manner of use of meters are subject to inspection in accordance with regulations and specifications established under the Electricity and Gas Inspection Act. The sealing and marking requirements are set forth in specifications established pursuant to section 18 of the Electricity and Gas Inspection Regulations. Installation and use requirements are set forth in specifications established pursuant to section 12 of the Regulations. Verification of conformity is required in addition to this approval for all metering devices excepting instrument transformers. Inquiries regarding inspection and verification should be addressed to the local inspection office of Industry Canada.

Original copy signed by:

Patrick J. Hardock, P.Eng.
Senior Engineer – Gas Measurement
Engineering and Laboratory Services Directorate

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ÉVALUÉ PAR**AG-0489 et Révision 3**
AG-0489 Révisions 1 et 2