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NOTICE OF APPROVAL – AVIS D'APPROBATION

METTLER ELECTRONIC TOP LOADING BALANCES PC400, PC400C, PC440, PC2000, PC4000, PC4400 and PCI6

BALANCES A PLATEAU EN TERE METTLER ELECTRONIC PC400, PC400C, PC440, PC2000, PC4000, PC4400 ET PCI6

Company: Fisher Scientific Ltd.,
112 Colonnade Road
Nepean, Ontario

Société: Fisher Scientific Ltd.,
112 Colonnade Road
Nepean, Ontario

Manufacturer: Mettler Instruments AG
CH-8606 Greifensee
Switzerland

Fabricant: Mettler Instruments AG
CH-8606 Greifensee
Suisse

Type of Device: Electronic balances that utilize electromagnetic force compensation to sense weight.

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Note: *Delta Range - this is a switch selective mode that increases the increment by a factor of 10 and reduces capacity by the same factor.

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All PC models have a Delta Display whereby the last decimal place is blanked out on initial application of weight but as the weight settles it re-appears. This function can be selected by a toggle switch located on the back of the housing. This

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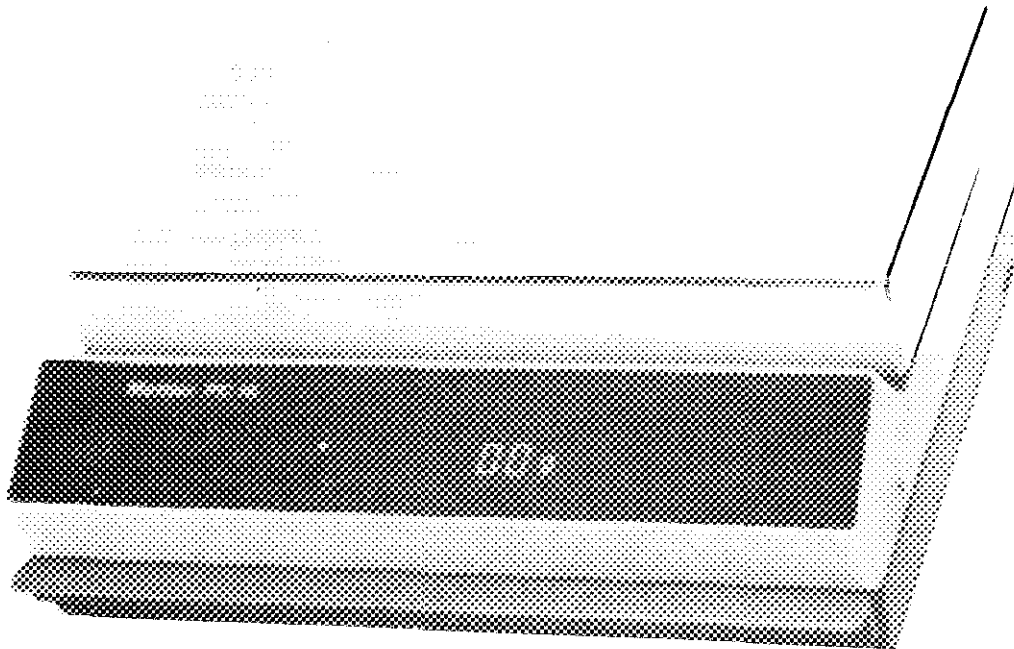
i.e.; on the PCI6I6 time for normal weight display is 0.75 sec. to 1.5 sec., with Delta Display ON 0.2 sec. to 0.4 sec.

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Approved Accessories: The following accessories are approved and shall be used with the models indicated:

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Application: For exclusive use in weighing precious metals and other commodities of comparable value.

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Description: These are top loading electronic balances. Weight is sensed using an electromagnetic force compensation principle. All models are fitted with levelling feet and a bulls-eye level. Some models must be operated with a draft shield as the weight increment and sensitivity are such that air currents will prevent the balance from having a true zero. These balances are fitted with + and - indication to the left of the weight display; when both are off, a zero display indicates that the zero position is accurate to $\pm \frac{1}{4}$ of an increment.

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There is only one operator control for these balances which controls, ON/OFF, zero and tare.

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To switch ON - press bar
To switch OFF - press bar
To zero - press bar. It operates full scale capacity

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To tare - press bar. It operates to full scale capacity.

Blanking for over capacity is indicated by a series of flat bars on the weight display. Sealing of the housing is by lead and wire seal through drilled head screws or Tamper proof paper seal.

Special Conditions:

1) As these balances have one display only, they must be labelled with instructions that state "THIS BALANCE MUST BE LOCATED SO THAT THE WEIGHT DISPLAY IS VISIBLE TO ALL PARTIES OF ANY TRADE TRANSACTION" (or words to that effect).

2) These balances must be installed on a solid, stable surface free from vibration and in an environment free from air currents.

3) These balances are exempted from the following requirements.

- a) Zero not to exceed 4% of capacity
- b) To identify "TARE" when a device is equipped to display only one weight value of a time
- c) To have levelling feet lockable

Conditions of Approval: Approval is granted under the Weights and Measures Act, S.C. 1970-71-72, chapter 36, and the Weights and Measures Regulations C.R.C.c., 1605 for use in Canada under the general conditions of the said Regulations, and under any special conditions listed above.

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