Ottawa **JAN 10 1983**

NOTICE OF APPROVAL – AVIS D'APPROBATION

Company: Ohaus Scale Corporation
29 Hanover Road
Florham Park, N.J.
USA 07932

Société: Ohaus Scale Corporation
29 Hanover Road
Florham Park, N.J.
USA 07932

Manufacturer: Ohaus Scale Corporation
29 Hanover Road
Florham Park, N.J.
USA 07932

Fabricant: Ohaus Scale Corporation
29 Hanover Road
Florham Park, N.J.
USA 07932

Type of Device: Dual range electronic
balance.

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Accessories: None

Accessoires: Aucun

Application: For exclusive use in
weighing precious metals and other
commodities of comparable value.

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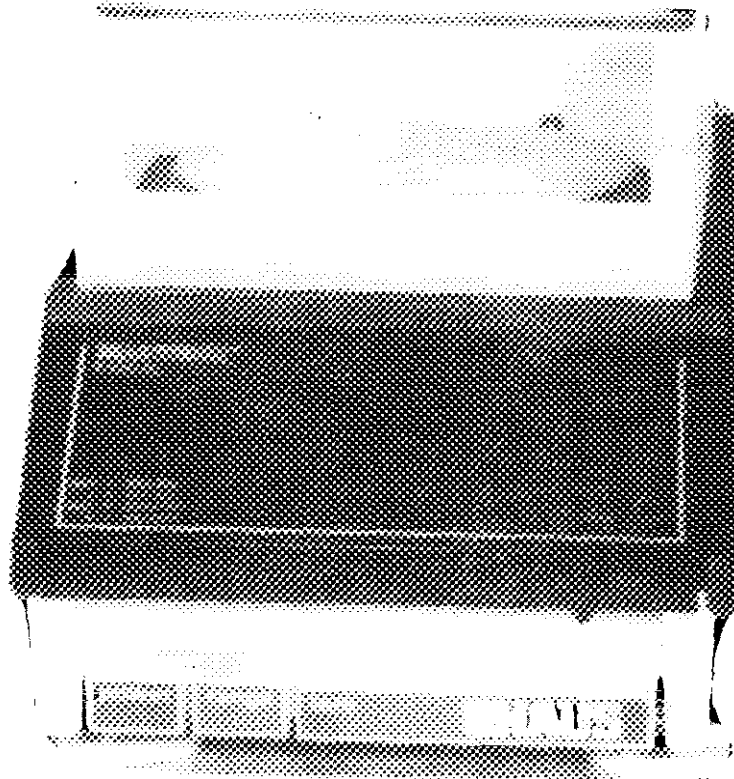
Description: These are dual range top
loading electronic balances. Weight is
sensed using an electromagnetic force
compensation principle. These models
must be operated with a draft shield as
the weight increment is such that air
currents will prevent the balance from
indicating a true zero.

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Display Switch: This switch is mounted
on the front of the balance directly
below the front window. This allows for
reading in grams or carats.

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- 2 -

Description: Continued

The following operator controls are located on the front of the scale:

- 1) "ON/OFF" power switch
- 2) "TARE" button has a dual purpose, it is used to tare off a weight on the platter, or it may be used for re-zeroing the scale.
- 3) DUAL RANGE OPERATION: the dual range offers both a high capacity and low capacity range with increased readability. To change ranges, start with the balance in the high capacity range. By depressing the range switch, the balance can be changed to the lower capacity range with increased readability. The changing of the range (high to low) will automatically tare the display which will produce a zero reading.

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Special Conditions:

- 1) These balances have one display only, they must be labelled with instructions that state "this balance must be located so that the weight display is visible to all parties of any trade transaction" (or words to that effect).
- 2) These balances are exempted from the requirements of SGM3/8.
- 3) These balances must be installed on a solid stable surface, free from vibration, and in an environment free from air current.

Condition of Approval: Approval is granted under the Weights and Measures Act, S.C. 1970-71-72, chapter 36, and the Weights and Measures Regulations, R.C.c., 1605 for use in Canada under the general conditions of the said Regulations, and under any special conditions listed above.

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W.R. Virtue

Chief
Legal Metrology Laboratories

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