SEP 24 1984

**NOTICE OF APPROVAL**

Company: Fisher Scientific Ltd.  
Division of Allied Canada Inc.  
112 Colonade Road  
NEPEAN, Ontario  
K2E 7L6

Manufacturer: Mettler Instruments A.G.  
CH-8606 Greifensee,  
Switzerland

Type of Device: Electronic balances that  
utilize electromagnetic force  
compensation to sense weight.

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C E 400

C E 150

Accessories: Compatible accessories that  
meet the requirements and perform in  
accordance with the Weights and Measures  
Act, Regulations and Specifications may  
be used.

Description: These are top loading  
electronic balances. Weight is sensed  
using an electromagnetic force  
compensation principle. All models are  
fitted with levelling feet and a bullseye  
level. Some models must be operated with  
a draft shield as the weight increment  
and sensitivity are such that air  
currents will prevent the balance from  
having a true zero. These balances are  
fitted with + and - indication to the  
left of the weight display; when both are  
off, a zero display indicates that the  
zero position is accurate  $\pm \frac{1}{4}$  of an  
increment.

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Société: Fisher Scientific Ltd.  
Division of Allied Canada Inc.  
112 Colonade Road,  
NEPEAN, Ontario  
K2E 7L6

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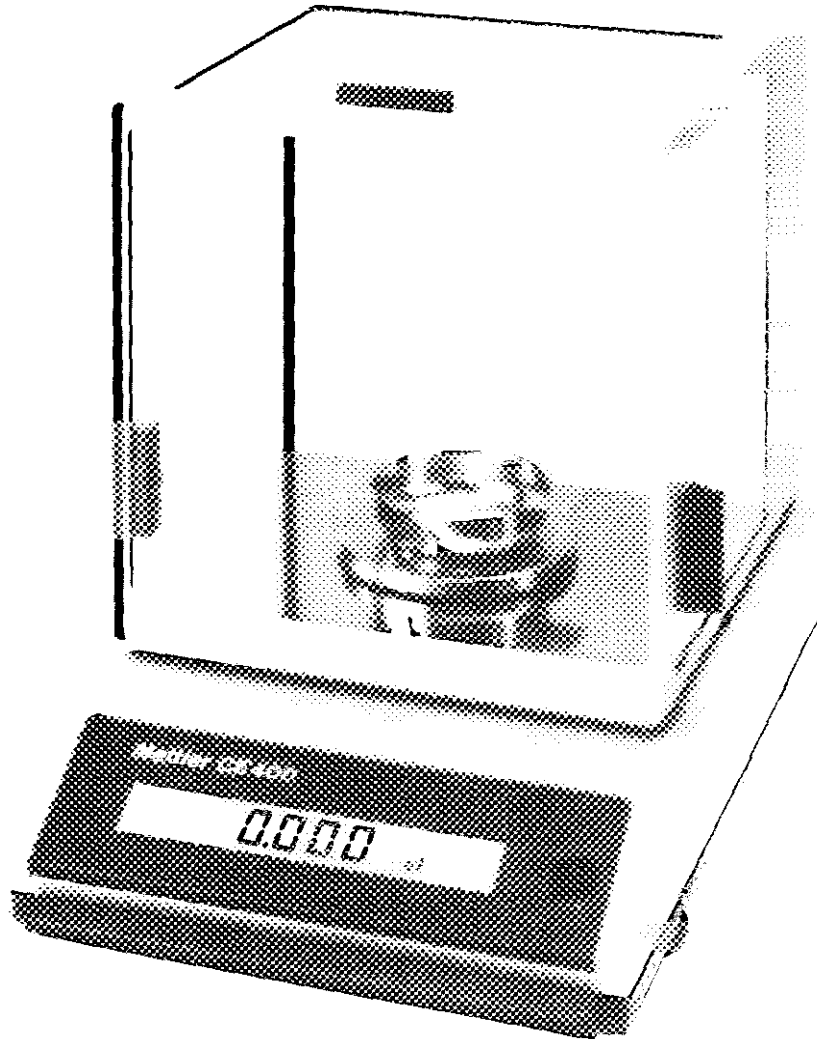
Description: Continued

There is only one operator control for these balances which controls, ON/OFF, zero and tare.

To switch ON - press bar

To switch OFF - lift bar

To zero - press bar.

Zero operates to full scale capacity.

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Description: Continued

To tare - press bar.  
Tare operates to full scale capacity.

Blanking for over capacity is indicated by a series of flat bars on the weight display. Sealing of the housing is by lead and wire seal through drilled head screws or tamper proof paper seal.

Special Conditions:

- 1) As these balances have one display only, they must be labelled with instructions that state "THIS BALANCE MUST BE LOCATED SO THAT THE WEIGHT DISPLAY IS VISIBLE TO ALL PARTIES OF ANY TRADE TRANSACTION" (or words to that effect).
- 2) These balances must be installed on a solid, stable surface free from vibration and in an environment free from air currents.
- 3) These balances are for exclusive use in weighing precious metals and other commodities of comparable value.
- 4) These balances are exempted from the following requirements:
  - a) Zero not to exceed 4% of capacity.
  - b) To identify "TARE" when a device is equipped to display only one weight value at a time.
  - c) To have levelling feet lockable.

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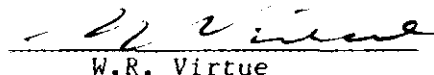
Conditions of Approval: Approval is granted under the Weights and Measures Act, S.C. 1970-71-72, chapter 36, and the Weights and Measures Regulations C.R.C.c., 1605 for use in Canada under the general conditions of the said Regulations, and under any special conditions listed above.

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W.R. Virtue

Chief  
Legal Metrology Laboratories

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