Consumer and
Corporate Affairs Canada

Legal Metrology

Consommation
et Corporations Canada

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APPROVAL No. - N° D'APPROBATION

S.WA-3212 Rev. 1

MS

AUG 21 1985

NOTICE OF APPROVAL

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Issued by statutory authority of
the Minister of Consumer and Corporate
Affairs under application by:

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NOTE: This approval applies only to
devices, the design, composition,
construction and performance of which
are, in every material respect,
identical to that described in the
information submitted and are typified
by the sample(s) submitted by the
applicant for evaluation for approval
in accordance with sections 14 and 15
of the Weights and Measures
Regulations. The following is a
summary of salient features only.

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SUMMARY DESCRIPTION:

These are top loading electronic balances. Weight is sensed using an electromagnetic force compensation principle. All models are fitted with lockable levelling feet and a bullseye level. These balances are fitted with + and - indication to the left of the weight display; when both are off, a zero display indicates that the zero position is accurate to $\pm \frac{1}{4}$ of an increment.

There is only one operator control for these balances which controls, ON/OFF, zero/tare and calibrate.

To switch ON - press bar

To switch OFF - lift bar

To zero - press bar. It operates to full scale capacity.

To tare - press bar. It operates to full scale capacity.

Blanking for over capacity is indicated by a series of flat bars on the weight display.

As these balances have one display only, they must be labelled with instructions that state "THIS BALANCE MUST BE LOCATED SO THAT THE WEIGHT DISPLAY IS VISIBLE TO ALL PARTIES OF ANY TRADE TRANSACTION" (or words to that effect).

These balances must be installed on a solid, stable surface free from vibration and in an environment free from air currents.

Zero is allowed to exceed 4% of capacity.

Tare is allowed without identification and visual confirmation that the tare mode is operating.

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SUMMARY DESCRIPTION: Continued

Note: *Delta Range - this is a mode that increases the increment by a factor of 10 and reduces capacity.

PE models with a Delta Display whereby the last decimal place is blanked out on initial application of weight but as the weight settles it re-appears. This function speeds up the weighing cycle.

Application - For exclusive use in weighing precious metals and other commodities of comparable value.

This Notice of Approval supersedes Notice of Approval S.WA-3212 dated October 15, 1984 and its addendum S.WA-3212 Add. 1 dated May 21, 1985.

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APPROVAL:

The design, composition, construction and performance of the device type(s) identified herein having been evaluated in accordance with regulations and specifications relating thereto, established under the Weights and Measures Act, approval is hereby granted pursuant to subsection 3(1) of the said Act.

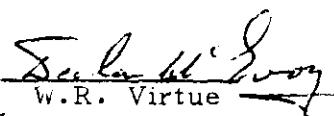
The marking, installation, use and manner of use in trade of devices are subject to inspection in accordance with regulations and specifications relating thereto, established under the Weights and Measures Act, and certification of conformity is required in addition to this approval. Inquiries regarding inspection and certification of conformity should be addressed to the local inspection office of Consumer and Corporate Affairs Canada. Requirements relating to marking are set forth in sections 18 to 26 of the Weights and Measures Regulations. Requirements relating to installation, use and manner of use are set forth in Part V and in specifications established pursuant to section 27 of the said Regulations.

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Chief
Legal Metrology Laboratories

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