Consumer and  
Corporate Affairs Canada

Consommation  
et Corporations Canada

Legal Metrology

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APPROVAL No. - N° D'APPROBATION  
S.WA-4049 Rev. 2

SEP 12 1988

**NOTICE OF APPROVAL**

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Issued by statutory authority of the  
Minister of Consumer and Corporate  
Affairs under application by:

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Fisher Scientific  
Division of Allied Canada Inc.  
112 Colonnade Road  
Nepean, Ontario  
K2E 7L6

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PM 34 Delta Range/Échelle Delta

0-4 kg x 0.1 g  
4-30 kg x 1 g

**NOTE:** This approval applies only to  
devices, the design, composition,  
construction and performance of which  
are, in every material respect,  
identical to that described in the  
information submitted; and are typified  
by the sample(s) submitted by the  
applicant for evaluation for approval  
in accordance with sections 14 and 15  
of the Weights and Measures  
Regulations. The following is a  
summary of principal features only.

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## SUMMARY DESCRIPTION:

These approved devices are electronic bench scales. Weight is sensed using an electromagnetic force compensation principle. They are fitted with lockable levelling feet and a bullseye level.

These devices are equipped with a single operator control which performs the following functions:

To switch ON - press bar.

To switch OFF - lift bar.

To Zero - press bar.

A graphic display "Delta Trac" on the right side of the weight display gives an idea of the used and unused portions of the weighing range.

Blanking for over capacity is indicated by a series of flat bars on the weight display.

These models incorporate a Delta Display whereby the last decimal place is blanked out on initial application of weight but as the weight settles it re-appears. This function speeds up the weighing cycle.

Delta Range - This is a mode that decreases the increment size by a factor of 10 and reduces capacity.

The Tare mode is allowed only when these devices are used in weighing precious metals and other commodities of comparable value, or determination of dockage in the grain industry. In this case, they are exempt from the requirements of sections 8, 9.2 and 9.6 of SGM3 and must be installed on a solid, stable surface free from vibration and in an environment free from air currents.

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## SUMMARY DESCRIPTION: Continued

If the installation does not permit customers to see the primary indicator, a remote indicator model GM48-2 or GM49-2 can be interfaced to the scale.

## APPROVAL:

The design, composition, construction and performance of the device type(s) identified herein have been evaluated in accordance with regulations and specifications established under the Weights and Measures Act. Approval is granted accordingly, pursuant to subsection 3(1) of the said Act.

The marking, installation, use and manner of use in trade of devices are subject to inspection in accordance with regulations and specifications established under the Weights and Measures Act. Verification of conformity is required in addition to this approval. Requirements relating to marking are set forth in sections 18 to 26 of the Weights and Measures Regulations. Requirements relating to installation, use and manner of use are set forth in Part V and in specifications established pursuant to section 27 of the said Regulations. Inquiries regarding inspection and certification of conformity should be addressed to the local inspection office of Consumer and Corporate Affairs Canada.

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W.R. Virtue

SEP 12 1988

Chief  
Legal Metrology Laboratories

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