MAR 3 0 1989

**NOTICE OF APPROVAL**

Issued by statutory authority of the Minister of Consumer and Corporate Affairs Canada for (category of device):

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**MODEL(S) / MODÈLE(S):**

Series/Série I...-\*K2  
Series/Série I...-XK2

**NOTE:** This approval applies only to devices, the design, composition, construction and performance of which are, in every material respect, identical to that described in the material submitted, and that are typified by samples submitted by the applicant for evaluation for approval in accordance with sections 14 and 15 of the Weights and Measures Regulations. The following is a summary of principal features only.

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**SUMMARY DESCRIPTION:**

The approved device is a top loading electro-mechanical balance. Weight is sensed using an electro-magnetic force compensation principle. The balance is fitted with levelling feet and a bullseye level. The levelling feet are fitted with a clamped-in insert of antisliding material at the top to prevent delevelling.

The LCD of 7 digits plus special characters provides a (+) or (-) indication to the left of the displayed weight value. When neither sign is lit, a zero display indicates that the zero position is accurate to  $\pm 1/4$  of an increment.

The operating program for this device is coded into an EEPROM. This internal switch which allows operator re programming of several parameters must be sealed to meet the requirements of SGM3/10. The keyboard "CAL" switch for external calibration must be disabled.

The device is equipped with operator controls which perform the following functions:

- ON/OFF; switches the display ON or OFF.
- CAL; to calibrate the balance by an internal motorized calibration device with a built-in weight piece. An internal switch locks this system for external calibration.
- TARE "T"; returns weight of empty container to zero balance. This key may also be used to switch display ON.

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Covered holes provide access to the calibration lock, linearity adjustment, and temperature resistor adjustment, and must be sealed.

This device is exempt from the requirements of sections 8, 9.2 and 9.6 of SGM3 when used as a pre-pack or check weigh scale.

Model Number Coding -  
Examples I...-\*K2; I...-XK2  
Where

- I = "Industry" model series
- ... = Capacity of balance (100g to 12100g)
- \*K2 = constructed for use in non-hazardous environments
- XK2 = constructed for use in hazardous environments.

#### APPROVAL:

The design, composition, construction and performance of the device type(s) identified herein have been evaluated in accordance with regulations and specifications established under the Weights and Measures Act. Approval is hereby granted accordingly pursuant to subsection 3(1) of the said Act.

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The marking, installation, use and manner of use in trade of devices are subject to inspection in accordance with regulations and specifications established under the Weights and Measures Act. Requirements relating to marking are set forth in sections 18 to 26 of the Weights and Measures Regulations. Installation and use requirements are set forth in Part V and in specifications established pursuant to section 27 of the said Regulations. A verification of conformity is required in addition to this approval. Inquiries regarding inspection and verification should be addressed to the local inspection office of Consumer and Corporate Affairs Canada.

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W.R. Virtue

Chief,  
Legal Metrology Laboratories

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