Consumer and
Corporate Affairs Canada
Legal Metrology

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MAY - 4 1992
MAI

NOTICE OF APPROVAL

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Issued by statutory authority of the Minister
of Consumer and Corporate Affairs Canada
for (category of device):

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NOTE: This approval applies only to
devices, the design, composition,
construction and performance of which are,
in every material respect, identical to that
described in the material submitted, and
that are typified by samples submitted by
the applicant for evaluation for approval in
accordance with sections 14 and 15 of the
Weights and Measures Regulations. The
following is a summary of principal features
only.

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Canada

SUMMARY DESCRIPTION:

The approved devices are top loading electronic balances. Weight is sensed using an electromagnetic force compensation principle. All models are fitted with lockable levelling feet and a bullseye level.

The devices are equipped with:

- Three single function keys.
 - OFF
 - PRINT (date output)
 - F (units, %, pcs)
- One multi function key
 - ON
 - RE-ZERO
 - MENU

Scales whose model numbers are suffixed with a "K" incorporate an internal calibration weight; pressing the re-zero key for a few seconds initiates a calibration sequence whereby the device autocalibrates. Calibration of other balances is effected similarly except calibration weights are added manually. With these scales, tare and re-zero are synonymous and both operate to full capacity.

The autocalibrate feature is not sealable; however, a change of greater than five graduations during autocalibration will abort the process.

The "Delta Trac" display is a circular dynamic graphic indicator that illuminates up to 60 radial segments in proportion to the load on the scale.

With the model PM 34-K Delta Range, pressing the re-zero bar with an applied load whose weight is between 0 and 32 000 g effects a 10-fold increase in readability within a window of $\pm 4\ 000$ g. The device will blank at 32 000 g.

With all scales, the last decimal place is suppressed during weighing, and will display only when the balance is in a stable condition.

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SUMMARY DESCRIPTION: Cont'd

The devices are for use in weighing precious metals and other commodities of comparable value, or determination of dockage in the grain industry, or for determining the net content of packages that are not weighed in the presence of the purchaser at the time of sale.

The devices are for use over a temperature range of +10°C to +30°C.

The units of weight are selectable from: ounce, pound, troy ounce, grain, carat, gram and kilogram.

The counting capability is not subject to Weights and Measures approval or inspection, but may be used in trade.

The devices are exempt from the requirements of sections 8, 9.2 and 9.6 of SGM3.

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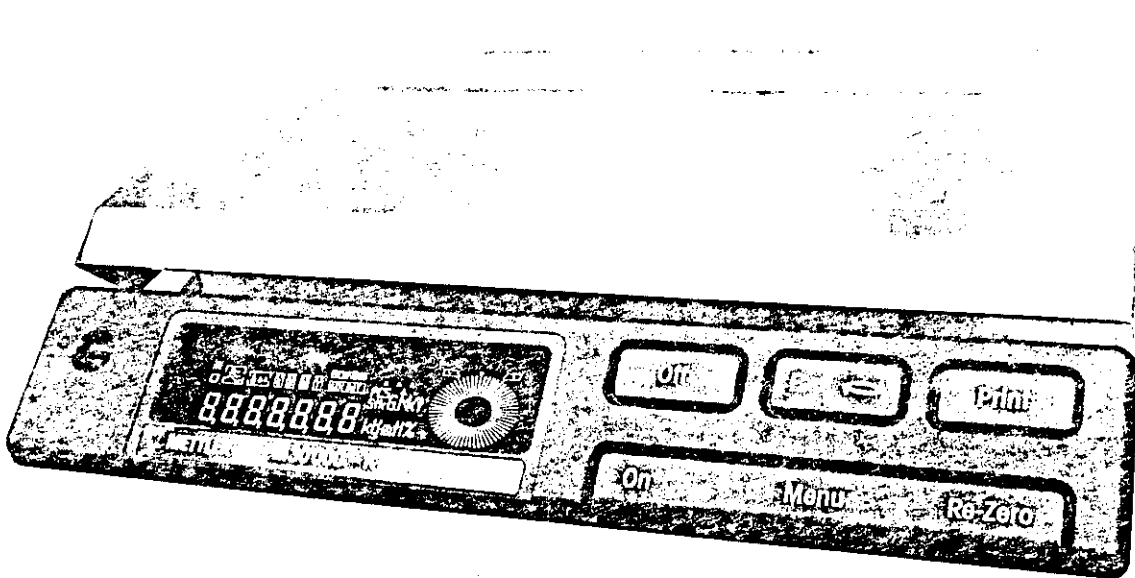
METTLER MODEL/Modèle PM 30000-K

APPROVAL:

The design, composition, construction and performance of the device type(s) identified herein have been evaluated in accordance with regulations and specifications established under the Weights and Measures Act. Approval is hereby granted accordingly pursuant to subsection 3(1) of the said Act.

The marking, installation, use and manner of use in trade of devices are subject to inspection in accordance with regulations and specifications established under the Weights and Measures Act. Requirements relating to marking are set forth in sections 18 to 26 of the Weights and Measures Regulations. Installation and use requirements are set forth in Part V and in specifications established pursuant to section 27 of the said Regulations. A verification of conformity is required in addition to this approval. Inquiries regarding inspection and verification should be addressed to the local inspection office of Consumer and Corporate Affairs Canada.

W.R. Virtue

Chief,
Legal Metrology Laboratories

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