SEP - 8 1995

NOTICE OF APPROVAL

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Issued by statutory authority of the Minister of Industry
for:

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RATING / CLASSEMENT

See "Summary Description" / Voir "Description Sommaire"

NOTE: This approval applies only to devices, the design, composition, construction and performance of which are, in every material respect, identical to that described in the material submitted, and that are typified by samples submitted by the applicant for evaluation for approval in accordance with sections 14 and 15 of the Weights and Measures Regulations. The following is a summary of the principal features only.

SUMMARY DESCRIPTION:

CATEGORY

The device is an electronic top loading bench scale featuring a liquid crystal display (LCD), three function keys and a stainless steel platter.

DESCRIPTION

A single, centrally located weight sensor, working on the electromagnetic force compensation principle, is encased within a plastic housing which is itself supported by one fixed and three lockable and adjustable feet. The device is powered by 12 VDC from a 110 VAC transformer, either directly or via an optional model PP-B10 rechargeable battery accessory.

DISPLAY

The LCD features an 8-digit, 7-segment indicator for weight display and alphabetic messages as well as icons and annunciators for, among others, motion and unit of weight.

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SUMMARY DESCRIPTION: Cont'd**FUNCTION KEYS**

Three function keys sublabelled "OFF", "F" and "Cal/Menu" provide the following functions;

- "OFF" switches the device on and off, effects platter tare/zero and aborts calibration and menu selection when in the Cal/Menu mode;
- "F", while 'Menu' is active, cycles among menu choices and
- "Cal/Menu" accesses, selects and stores values when in the calibrate/menu mode.

CALIBRATION

- a) If the device is equipped with a built-in calibration weight, autocalibration can be effected by keyboard entry. This calibration function is not sealable; however, a change of greater than five "d" during calibration will abort the process.
- b) For devices not so equipped, the calibration routine is accessed via a sealable device setup mode (refer to SEALING). The user is prompted to unload and then load a specified test weight at which point the device autocalibrates.

SEALING

Device setup is enabled:

- for the PB series, by removing, rotating 180° and reinserting an access button located on the right side of the device and
- for the SB series, by depressing a calibration switch located beneath the platter.

Proper sealing requires tamper proof seals to be placed over the access button (PB series)/calibration switch (SB series) and over the seam at the back of the device, sealing the two housing halves. The design is exempt from providing ready access to all other components or adjustments without breaking a seal as specified in SGM 3/10.

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SUMMARY DESCRIPTION: Cont'd**DELTARANGE**

For devices equipped with the DeltaRange feature, depressing the zero/tare function key with a load on the platter effects a 10-fold increase in readability within a reduced range (see Summary Chart).

COUNTING CAPABILITY

The ability to count, although not subject to Weights and Measures approval and initial inspection, may be used in trade.

LEVEL CONDITION

Device level is ensured by means of three lockable and adjustable levelling feet and a bull's eye level.

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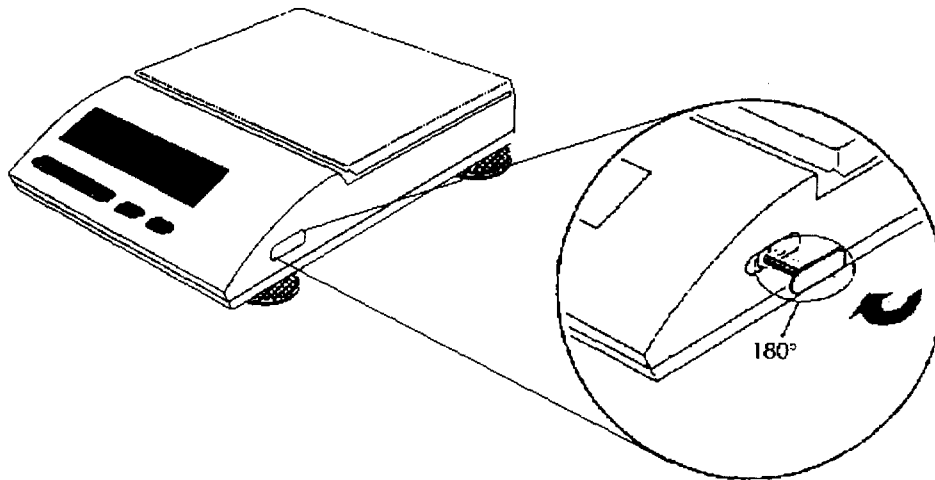
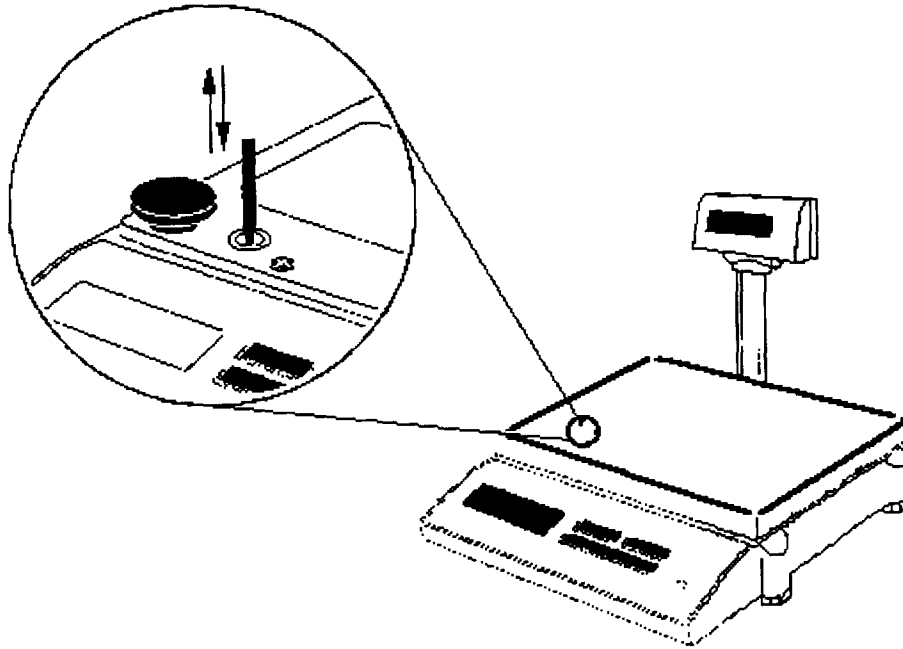
Series / série SB

USE RESTRICTION

The device is for use exclusively in weighing precious metals and other commodities of comparable value, or in the grain industry, or the determination of the net content of packages that are not weighed in the presence of the purchaser at the time of sale.

TEMPERATURE RANGE

The device is approved for use within a temperature range of +10 °C to +30 °C.

UNITS OF WEIGHT

The device is approved for use with the following units: kilograms, milligrams, pounds, ounces, troy ounces, grains and carat. Depending on the model, any two units may be chosen while in the setup mode and subsequently selected by the user via the "F" key.

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SUMMARY DESCRIPTION: Cont'd**VERIFICATION INTERVAL**

The value of the scale division (displayed), "d", which is the difference between two consecutive indicated values, is differentiated from the value of the verification scale division, "e", by square brackets about "d" within the display. For approval and inspection purposes, the limits of error applicable to the device are based on the value of "e".

EXEMPTION

Zero and tare being synonymous with these devices, they are thus exempt from the requirements of sections 8 and 9.6 of SGM 3.

REVISION

The purpose of Revision 1 was to add the GB Series and corrections to Pan dimensions.

The purpose of Revision 2 is to correct the pan dimensions in the "Model Summary Chart" on page 7.

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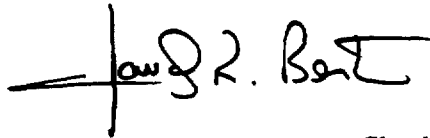
APPROVAL:

The design, composition, construction and performance of the device type(s) identified herein have been evaluated in accordance with regulations and specifications established under the Weights and Measures Act. Approval is hereby granted accordingly pursuant to subsection 3(1) of the said Act.

The marking, installation and manner of use of trade devices are subject to inspection in accordance with regulations and specifications established under the Weights and Measures Act. Requirements relating to marking are set forth in sections 18 to 26 of the Weights and Measures Regulations. Installation and use requirements are set forth in Part V and in specifications established pursuant to section 27 of the said Regulations. A verification of conformity is required in addition to this approval. Inquiries regarding inspection and verification should be addressed to the local inspection office of Industry Canada.

Claude R. Bertrand, P.Eng.
A/Manager
Approval Services Laboratory

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