**NOTICE OF APPROVAL**

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Issued by statutory authority of the Minister of Industry for:

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**NOTE:** This approval applies only to devices, the design, composition, construction and performance of which are, in every material respect, identical to that described in the material submitted, and that are typified by samples submitted by the applicant for evaluation for approval in accordance with sections 14 and 15 of the Weights and Measures Regulations. The following is a summary of the principal features only.

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## SUMMARY DESCRIPTION:

### CATEGORY

The device is an electronic portable bench scale (balance).

### DESCRIPTION

The approved device is a self-contained electronic balance, that is fitted with a single centrally located weight sensor, working on the electromagnetic force compensation principle. The weight sensor supports a stainless steel load receiving element. The device components are encased within an aluminum housing which is supported by three (3) feet, one fixed (or stationary) in the front center of the device and two (2) adjustable in each rear corner. The device can be fitted with a draft shield and is powered by a 12 VDC from a 100 VAC external transformer.

### DISPLAY

The device features a vacuum fluorescent indicator for weight display, alphabetic messages, delta range graphic, weighing ranges, delta range for finer readability and annunciators for among others, motion, cal, load, done, and abort.

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## FUNCTION KEYS

The device is fitted with the following keys:

“ON/OFF”, “TARE”, “e”, “←0← cal”, “F” and “Menu”.

Note that the zero and tare key functions (*O/T*) are synonymous with these devices.

## CALIBRATION

The device can be calibrated automatically with a built-in calibrated weight. In this mode, the balance will perform an automatic calibration sequence when needed. During this automatic calibration sequence the display will show the words “CAL”. When this sequence is complete, the device will return to normal operation.

## SEALING

The scale can be sealed with a wire security seal threaded through a hole in the top and bottom halves of the rear of the enclosure preventing the housing from being separated. It is also necessary to place a self-destructive seal over one of the screws under the sub-platter to prevent access to calibration and configuration parameters.

## SERIAL CONNECTIONS

Standard design RS232C

## TEMPERATURE RANGE

The device is approved for use within a temperature range of 5°C to 40°C.

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Suffix “DR”: Delta-Range: The displayed division “d” changes at the lower capacity listed for the model. The verification division “e” remains constant.

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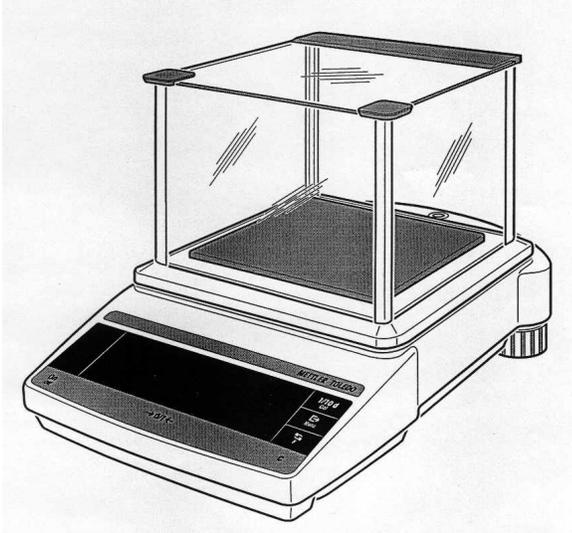
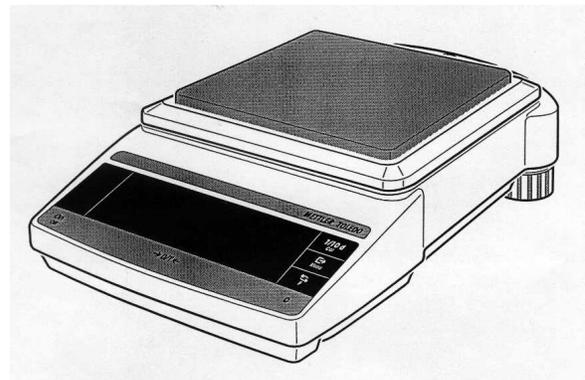
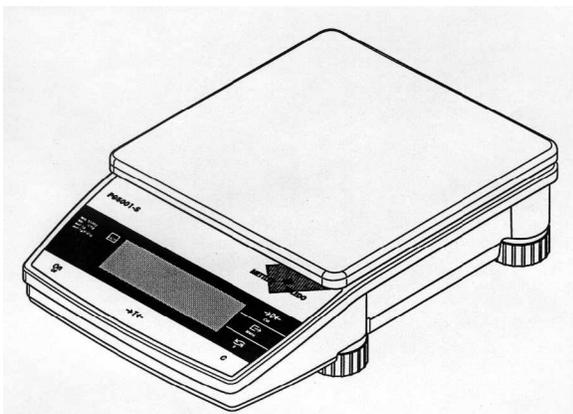
240 mm x/sur 240 mm Platter / plateau

**APPROVAL:**

The design, composition, construction and performance of the device type(s) identified herein have been evaluated in accordance with regulations and specifications established under the Weights and Measures Act. Approval is hereby granted accordingly pursuant to subsection 3(1) of the said Act.

The marking, installation and manner of use of trade devices are subject to inspection in accordance with regulations and specifications established under the Weights and Measures Act. Requirements relating to marking are set forth in sections 49 to 54 of the Specifications Relating to Non-Automatic Weighing Devices. Installation and use requirements are set forth in sections 55 to 67 of the Specifications Relating to Non-Automatic Weighing Devices. A verification of conformity is required in addition to this approval. Inquiries regarding inspection and verification should be addressed to the local inspection office of Industry Canada.

Original copy signed by:

René Magnan, P. Eng  
Director  
Approval Services Laboratory

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