NOTICE OF APPROVAL

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Issued by statutory authority of the Minister of Industry
 for:

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APPLICANT

REQUÉRANT

McKesson Automated Prescription Systems
 4333 Shreveport Hwy
 Pineville, Louisiana, 71360
 USA

MANUFACTURER

FABRICANT

McKesson Automated Prescription Systems
 4333 Shreveport Hwy
 Pineville, Louisiana, 71360
 USA

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RATING/ CLASSEMENT

BU 2010

Max

d

e_{min}

n_{max}

620 g

0.001 g

0.01 g

62 000

Accuracy Class / Categorie de precision: II

NOTE: This approval applies only to devices, the design, composition, construction and performance of which are, in every material respect, identical to that described in the material submitted, and that are typified by samples submitted by the applicant for evaluation for approval in accordance with sections 14 and 15 of the Weights and Measures Regulations. The following is a summary of the principal features only.

SUMMARY DESCRIPTION:

CATEGORY

The approved device is an electronic self-contained bench scale.

DESCRIPTION

The device is powered by an AC/DC adapter.

Weight is sensed by an electromagnetic force compensating load cell which is centrally located. The load cell is bolted to an aluminium base and supports a spider assembly with a stainless steel platter measuring 135 mm x 135 mm.

The capacity of the load cell is 620 g and is protected from overloading by means of overload stops. The device is levelled by means of two adjustable feet and a bull's eye level.

The device is housed in a cast aluminium casing and utilises a digital vacuum fluorescent display.

Annunciators are: LB/G (gross), PT (tare), NET, $\div 0^2$ (center of zero), \mathbf{y} (wt is below zero) and t (weight exceeds zero) and others. The device has a counting feature which is not legal for trade.

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CONTROL KEYS

The operator controls are performed via a 10 button keypad and permit the following choices among others:

ON/OFF RE-ZERO	used to turn the scale on or off or to rezero the scale
MENU/esc	used to exit the menu or to enter application and configuration menu
⇌ ÷	used to change from menu path into sub-paths and back or to select from respective display
€ •	used to move up/down within the main and sub-paths or to select from respective display
T/5	used to select parameters, stores the changed parameters or to tare the scale
T/ins	used to tare the scale or to place insert markers in text entry
MODE/Clr	used to delete input (in text entry), to change mode from Weighing, Manual Count and APS application
RxCheck/Start	used to reset to beginning of APS application (except in Host Mode).

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10°C to 30°C / 50°F to 86°F)

SEALING

Means of adjustment and configuration parameters can be sealed in two ways or a combination of using both methods:

1. Two drilled head screws, one that secures the top and bottom sections of the device and the other drilled screw that when inserted into the hole (located just below the screw that secures the top and bottom pieces) will not allow access to the calibration button.
2. With a screw already inserted into the calibration hole, a flat tapered head screw with a long shaft that would extend down past the calibration access screw, thus not allowing it to be backed out. This flat, tapered screw when fully inserted can be sealed with a paper seal covering the head of the screw and onto the metal sides surrounding it.

EVALUATED BY

Ken Chin
Complex Approvals Examiner
Tel: (613) 954-2481
Fax: (613) 952-1754

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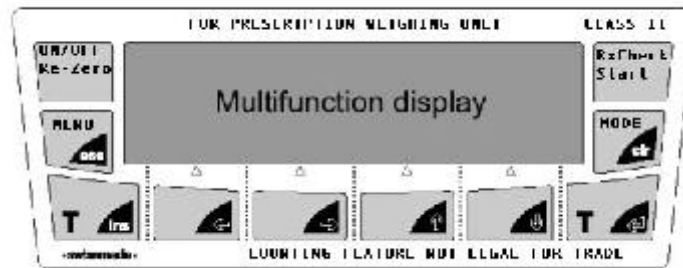
APPROVAL:

The design, composition, construction and performance of the device type(s) identified herein have been evaluated in accordance with regulations and specifications established under the Weights and Measures Act. Approval is hereby granted accordingly pursuant to subsection 3(1) of the said Act.

The marking, installation and manner of use of trade devices are subject to inspection in accordance with regulations and specifications established under the Weights and Measures Act. Requirements relating to marking are set forth in sections 49 to 54 of the Specifications Relating to Non-Automatic Weighing Devices. Installation and use requirements are set forth in sections 55 to 67 of the Specifications Relating to Non-Automatic Weighing Devices. A verification of conformity is required in addition to this approval. Inquiries regarding inspection and verification should be addressed to the local inspection office of Industry Canada.

Original copy signed by:

René Magnan, P. Eng
Director
Approval Services Laboratory

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